VersaTube (VT502)

Instructions for Use

1. Lift tab. Push down on lid if tab is stuck.
2. Open lid.
3. Fill as required.
4. Close lid to required venting angle.
5. Process sample.
6. If lid needs to be opened during processing, tap top of lid to remove excess condensed moisture before lifting lid.
7. To fully close lid for storage or centrifugation, use the palm of hand to firmly press lid completely flat and close tab.
8. To remove cap: twist or cut cap at hinge.

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STATEMENT OF ACCURACY - VOLUME GRADUATIONS FOR VT502 - VERSATUBE

THIS CERTIFIES THAT THE DISPOSABLE TUBES DEVELOPED FOR METALS DIGESTION HAVE BEEN TESTED BY AN INDEPENDENT LABORATORY IN THE USA TO DETERMINE THE ACCURACY OF VOLUME GRADUATIONS WHEN COMPARED TO CLASS A SPECIFICATIONS. TESTING IS BASED ON NIST SOP14 AND ASTM E1272. SAMPLES WERE FOUND TO BE ACCURATE TO +/- 0.25ML AT 50ML. TUBES WERE TESTED AT 20-23C AND DONE ON A "TO CONTAIN" BASIS.

METALS SCREENING STATEMENT FOR VT502 - VERSATUBE

A REPRESENTATIVE SAMPLE OF OF TUBES FROM A TYPICAL PRODUCTION LOT WAS TESTED FOLLOWING SOP 7040, REV 10 WHICH IS AMENABLE TO EPA METHOD 200.8 BY AN INDEPENDENT, ACCREDITED LAB IN THE USA. TESTS WERE PERFORMED BY ICPMS AND SAMPLES CHECKED FOR 66 ANALYTES. ALL TUBES WERE FOUND TO HAVE NON DETECTABLE LEVELS OF THE ANALYTES. IT IS IMPORTANT TO NOTE HOWEVER, THAT DEPENDING ON HOW THE TUBES ARE HANDLED, THE ENVIRONMENTAL CONDITIONS ETC. SUCH RESULTS MAY VARY - CUSTOMERS ARE ENCOURAGED TO RANDOMLY CHECK SAMPLES ON THEIR OWN TO ENSURE THEY MEET WITH THEIR INTERNAL REQUIREMENTS.

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